



# REFERENCE CARD for Charm® Chloramphenicol Test

This summary test procedure is not for use in regulatory testing. Refer to the current Operator's Manual for the complete procedure.

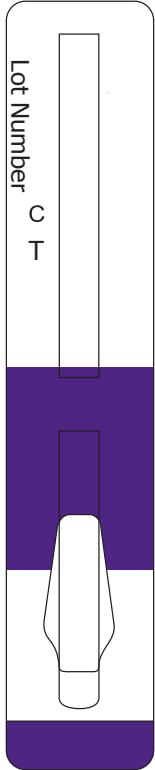


Place completed strip here.

## Invalid

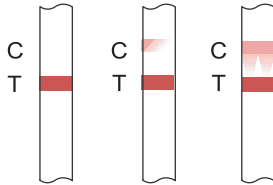


Re-test **Invalid** samples


## Positive

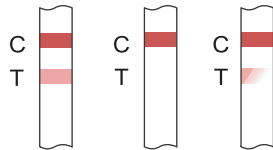

## Negative

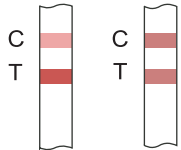

1

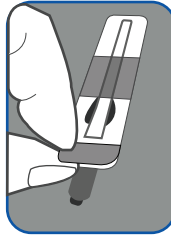


Verify incubator temperature is  $40 \pm 2$  °C.

Shake milk sample (or controls) and label test strip.

Place strip in incubator and use tab to peel tape back to "Peel to here" line.

## Peel Strip

2

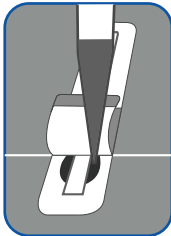

Slowly pipet 300  $\mu$ L sample (or controls) into sample compartment at ROSA® Incubator indicator line (as shown).

Reseal tape over sample compartment.

Close lid and latch.

Incubate for 8 minutes, but not more than 10 minutes.

## Add Raw Milk Sample

3



## Incubate

Remove tests from incubator.

Visually inspect C (control) line for even development.

Insert clean and visually valid test strip in ROSA Reader or Charm EZ® system in Read-only mode and read results.

T +1.978.687.9200 | F +1.978.687.9216 | E info@charm.com | www.charm.com



659 Andover Street, Lawrence, MA 01843-1032, USA

© 2016 Charm Sciences, Inc. Charm, Charm EZ and ROSA are registered trademarks of Charm Sciences, Inc.

MRK-258-003 Feb-2016