

mycotube

The 5 minute Aflatoxin field test you can use anywhere.

The **Mycotube Portable Aflatoxin Screening Test** is easy and fast:

- » No equipment, no solvent or electricity required
- » Detection levels between 10 to 100 ppb
- » Results in only 5 minutes

To get a clear and accurate test anywhere, you only need MycoTube, your sample, and water. Follow these six simple steps:

1



Put ground sample in MycoTube.

2




Add water.

3



Shake and settle.

4

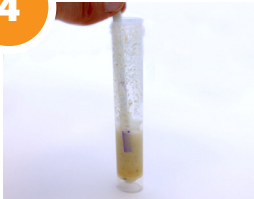

Add MycoTube strip.

5



Let sit for 5 minutes.

6



Read results.



About Charm

Charm is a leading manufacturer of food safety products. We create innovative solutions for mycotoxin, water quality, antibiotics, sanitation hygiene, and microbial detection. We offer practical and accurate solutions for use in field, plant, and laboratories.

Charm, Where the Science of Food Safety is a Way of Life.

Mycotube is the only rapid, qualitative visual assay to detect aflatoxin in maize and peanuts. Each MycoTube vial contains extraction powder and one lateral flow strip in dipstick format. MycoTube's cap doubles as a scoop for ground samples and water. Incubation is at room temperature. Designed to generate positives at 10 ppb.

For more information, or to order a free test kit, visit www.charm.com/FreeMycotube.

www.charm.com

T +1.978.687.9200 | F +1.978.687.9216 | E info@charm.com | 659 Andover Street, Lawrence, MA 01843
© 2018 Charm Sciences, Inc. Charm is a registered trademark of Charm Sciences, Inc. | MRK-1269-001 Oct-2018

