

mycotube

For Aflatoxin

The 5 minute Aflatoxin field test you can use anywhere.

The **Mycotube Portable Aflatoxin Screening Test** is easy and fast:

- » No equipment, solvent, or electricity required
- » Detection levels >10 ppb
- » Results in only 5 minutes

To get a clear and accurate test anywhere, you only need the MycoTube vial, your sample, and water. Follow these six simple steps:



1
Add ground sample into MycoTube vial.



2
Add water.




3
Shake and settle.



4
Insert MycoTube strip.



5
Let sit for 5 minutes.

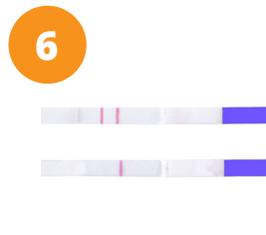

6
Read results.



About Charm

Charm is a leading manufacturer of food safety products. We create innovative solutions for mycotoxin, water quality, antibiotics, sanitation hygiene, and microbial detection. We offer practical and accurate solutions for use in field, plant, and laboratories.

Charm, Where the Science of Food Safety is a Way of Life.

The Mycotube for Aflatoxin test is the only rapid, qualitative assay to visually detect aflatoxin in maize and peanuts. Each MycoTube vial contains extraction powder and one lateral flow strip in dipstick format. The MycoTube cap doubles as a scoop for ground samples and water. The incubation is at room temperature and it is designed to generate positives above 10 ppb.

For more information, or to order a free test kit, visit www.charm.com/FreeMycotube.

www.charm.com

T +1.978.687.9200 | F +1.978.687.9216 | E info@charm.com | 659 Andover Street, Lawrence, MA 01843
© 2019 Charm Sciences, Inc. Charm is a registered trademark of Charm Sciences, Inc. | MRK-1269-001 Aug-2019

